

## **MICRONIC | SAMPLE HANDLING | CAPPING & DECAPPING**

### **UNIVO ELECTRIC CAPPER CP480**

**MAKE YOUR CAPPING EASY. MICRONIC OFFERS AN ELECTRIC CAPPER FOR THE SEALING OF A VARIETY OF TUBE SIZES WITH CAPMATS, CAPBANDS AND CAPCLUSTERS. THE ELECTRIC CAPPER SEALS AN ENTIRE RACK OF 96 TUBES WITHIN 7 SECONDS AND HAS A GUARANTEED EVENLY CAPPING RESULT.**



Built to meet the special individual requirements of our customers, the Univo Electric Capper can cap all Micronic plug style caps: capmats, capbands and capclusters made of EVA or TPE, on Micronic 96-format tubes and racks. The capper is also compatible with other brands of caps and tubes (please contact us for a list of compatible brands).

The Univo Electric Capper has a compact and robust design. For your convenience, the capper operates without any buttons. Simply pull the drawer out of the capper, place a rack filled with tubes on the drawer, and then place the capmat, capband or capcluster on the tubes. Push the drawer inside the capper and the capping will start automatically.

The controlled pressure system will press all the caps down evenly to ensure the integer capping of your tubes. Simply pull the drawer out and retrieve the rack. One capping cycle lasts approximately 7 seconds.

The Univo Electric Capper has the flexibility to cap any tube height up to 48mm (no adapters are required), and both full and partially-full racks of tubes. The pressure system of the capper is adjustable which ensures correct capping.

For more information:  
e-mail to [sales@micronic.com](mailto:sales@micronic.com)


**Micronic Europe B.V.**  
(Asia, Africa, Europe and Oceania)  
Phone: +31 (0)320 277 070


**Micronic America LLC**  
(North, Central and South America)  
Phone: +1 484 480 3372

Please visit our website: [www.Micronic.com](http://www.Micronic.com)

#### **INFORMATION**

- No adapters required
- Caps an entire rack of 96 tubes in approximately 7 seconds
- Indication light underneath when switched on
- Operates without any buttons
- Compatible with all Micronic tube sizes up to 1.40ml
- Also compatible with other brands of tubes and caps (please contact us for a list of compatible brands)
- Compact and robust design
- Controlled pressure system
- Works on full and partially-full racks

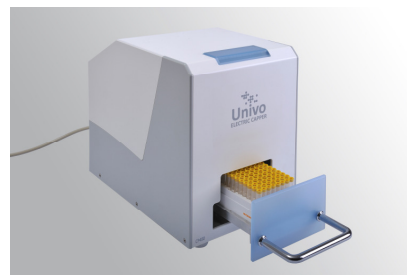

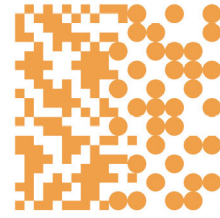

**MICRONIC**

## **MICRONIC | SAMPLE HANDLING | CAPPING & DECAPPING**

### **SPECIFICATIONS**

COMPATIBILITY	ALL SBS STANDARD 96-FORMAT RACKS
CYCLE TIME	7 SECONDS PER RACK
DIMENSIONS	38CM X 21CM X 29CM / 15INCH X 8.3INCH X 11.4 INCH
WEIGHT	14 KG / 31 LBS
VOLTAGE	115 V - 230 V



### **ORDERING INFORMATION**

#### **Univo Electric Capper CP480**

MP35020C Univo Electric Capper CP480 for all tubes sizes up to 1.40ml unit



Micronic has a broader product range than the products listed under ordering information. For our complete product portfolio, please send your e-mail to [sales@micronic.com](mailto:sales@micronic.com) or contact us by phone.



**MICRONIC**

**SAMPLE HANDLING | EQUIPMENT | M505.2 6.2**